



*Thin Film Measurement solution
Software, sensors, custom development
and integration*

TFCompanion Update to ver. 4.5

These instructions are only for the update from the previous versions of TFCompanion. If you already have ver. 4.5 of TFCompanion – updates to new builds are done as before: Help/Check Update and follow prompts.

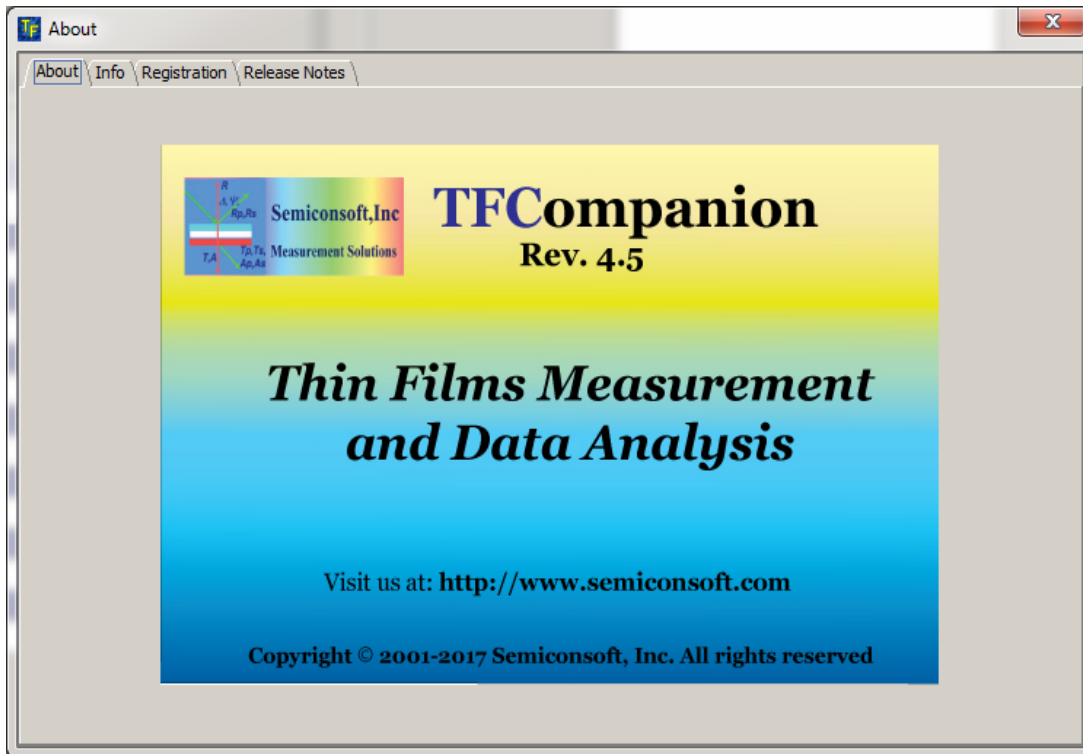


Fig. 1 Help/About from the main menu will show the current version of TFCompanion.

- I. Steps to update TFCompanion to ver. 4.5
 1. Install the new update (TFUpdater)
 - a). Most likely, you have a previous version of TFUpdater

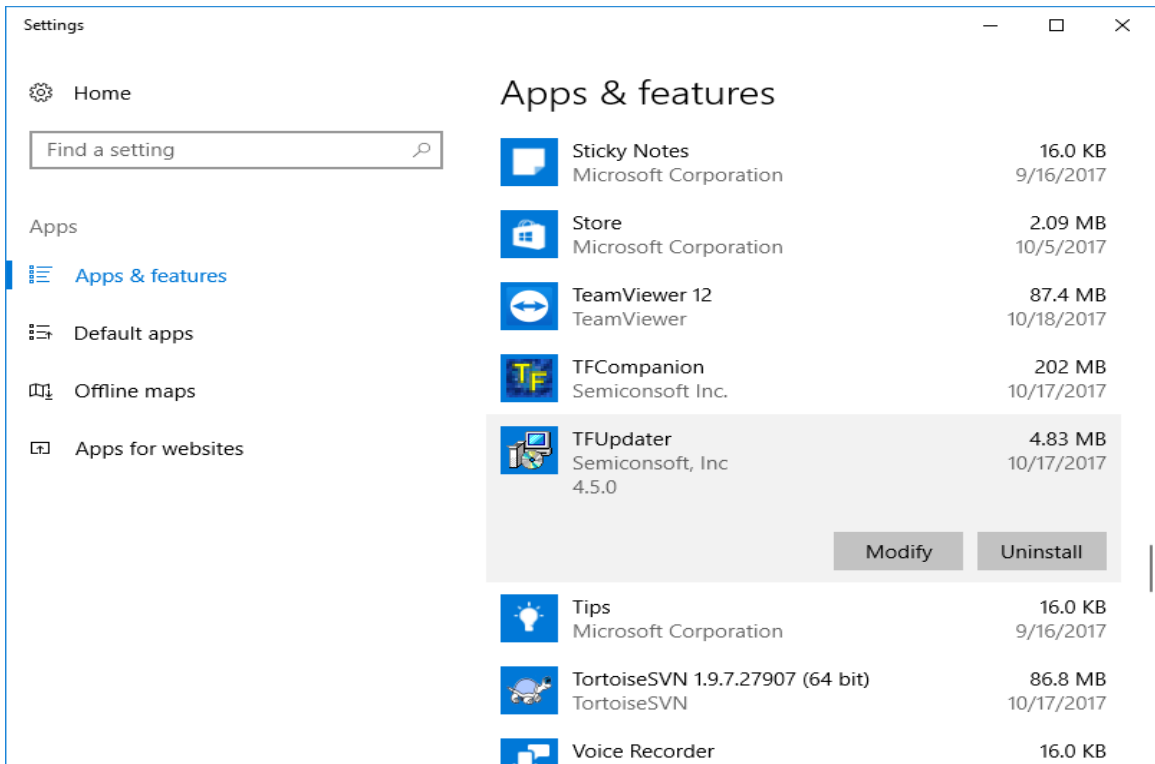


Fig. 2 Check the version of TFUpdater. If the version is < 4.5 – uninstall it

b). After uninstalling TFUpdater – delete C:/TFCompanion/updater directory

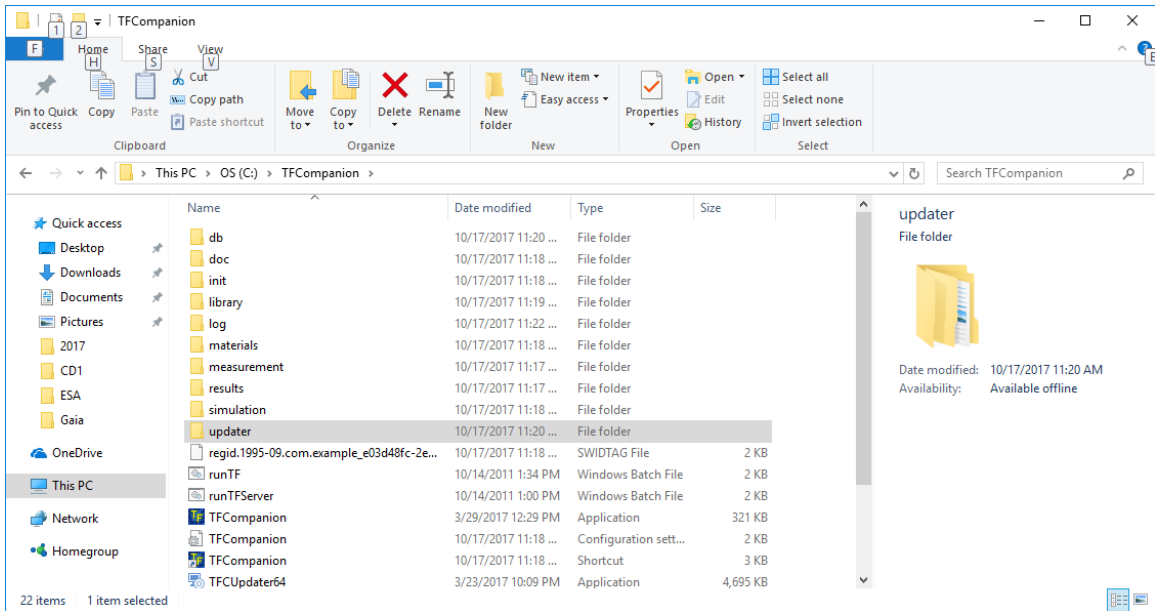


Fig. 3 C:/TFCompanion/updater directory need to be removed

c). Download and install new TFUpdater

64 bit version:

<http://www.semiconsoft.com/html/download/Installers/InstData/Windows/updater/TFUPdater64.exe>

32 bit version:

<http://www.semiconsoft.com/html/download/Installers/InstData/Windows/updater/TFUPdater.exe>

d). After new TFUpdater is installed – you are ready to update the software.
Start TFCompanion and use Help/Check update

2. If you are using Dynamic measurement plugin (Kinetic measurement). One of the libraries need to be updated manually.

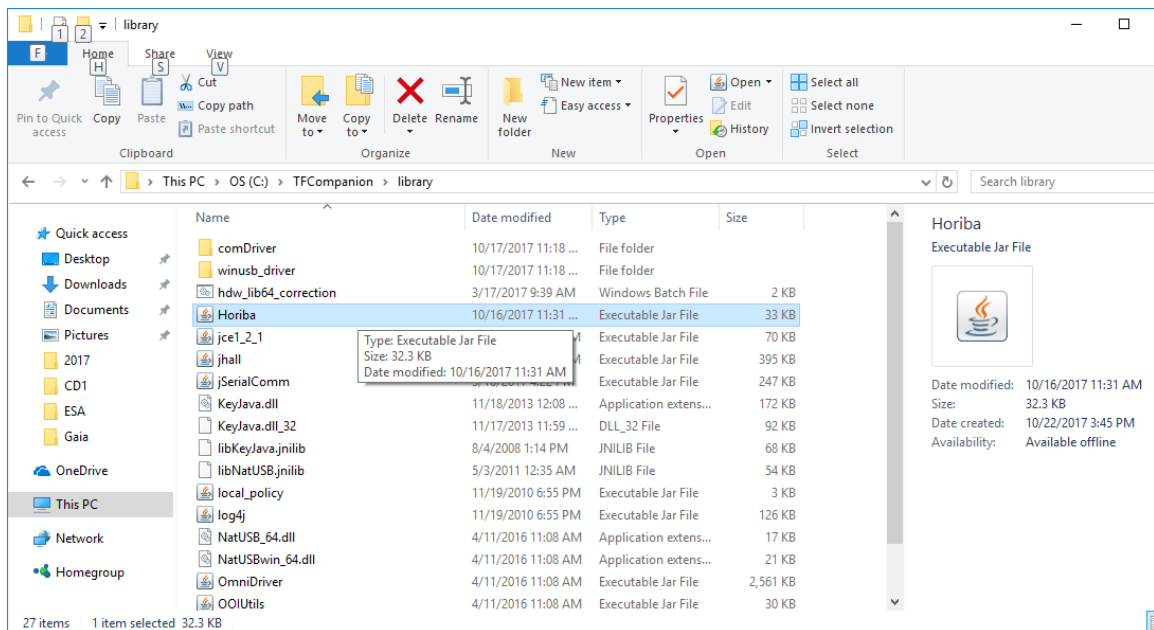


Fig. 4. Download file Horiba.jar: <http://www.semiconsoft.com/Exchange/Horiba.jar> and place it in the C:/TFCompanion/library directory